thermofisher.com

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☐ [Rn] Rat ⓘ

Enter target information keyword(s)
(Max 25 entries allowed)

Enter one or more keywords. Search multiple keywords by separating with a space, comma, or press Enter after each search keyword.

Supported Keywords: Assay IDs, NCBI Gene IDs, Gene Symbols, RefSeq Accession numbers, RefSeq Protein IDs.

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The screenshot shows the ThermoFisher siRNA search interface. At the top, it asks "What type of siRNA do you want ?" with four options: Silencer® Select siRNA (Recommended), Ambion® In Vivo siRNA, Silencer® siRNA, and Stealth™ RNAi siRNA. Below this, it asks "What species do you want to target ?" with three options: [Hs] Human, [Mm] Mouse, and [Rn] Rat. The third option is highlighted with a red box and a red circle with the number 1. Below the species selection, it asks "Enter target information keyword(s)" with a text input field and a "Search siRNA" button. The input field is highlighted with a red box and a red circle with the number 2. The "Search siRNA" button is highlighted with a red box and a red circle with the number 3. A red arrow points from the text box on the right to the input field.

What type of siRNA do you want ?

Recommended

☐ Silencer® Select siRNA [i](#) ☐ Ambion® In Vivo siRNA [i](#) ☐ Silencer® siRNA [i](#) ☐ Stealth™ RNAi siRNA [i](#)

What species do you want to target ?

☐ [Hs] Human  ☐ [Mm] Mouse  ☐ [Rn] Rat 

Enter target information keyword(s)
(Max 25 entries allowed)

Enter one or more keywords. Search multiple keywords by separating with a space, comma, or press
Enter after each search keyword.

Supported Keywords: Assay IDs, NCBI Gene IDs, Gene Symbols, RefSeq Accession numbers,
RefSeq Protein IDs.

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☐	Assay ID	Gene	Transcripts	Species	Transcript Type	Product Type	Specifications	Price (JPY)
☐	s8403 ① ★ Validated ②	SMAD4	6	Human	Coding ③	Silencer Select	Desalted 5 nmol ▼ ④	定価価格 Each CAT # 4390824 Estimated ship by: 4-5 business days Add to cart
View Details ▼ Genomic Map ▼ Validation Data ▼ Recommended Controls ▼ Recommended Products ▼								

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☐	Assay ID	Gene	Transcripts	Species	Transcript Type	Product Type	Specifications	Price (JPY)
☐	s8403	SMAD4	6	Human	Coding	Silencer Select	Desalted 5 nmol ▼	定価価格 Each CAT # 4390824 Estimated ship by: 4-5 business days
★ Validated								
View Details ▼ Genomic Map ▼ Validation Data ▼ Recommended Controls ▼ Recommended Products ▼								

Confirm Size & Purification

Please confirm your Size & Purification selections before adding to your cart

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☐ Assay ID	Gene	Transcripts	Species	Transcript Type	Product Type	Specifications	Price (JPY)
☐ s8403 ★ Validated	SMAD4	6	Human	Coding	Silencer Select	Desalted 5 nmol ▼	定価価格 Each CAT # 4390824 Estimated ship by: 4-5 business days
① View Details ▼ ② Genomic Map ▼ ③ Validation Data ▼ Recommended Controls ▼ Recommended Products ▼							

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Gene Symbol:	SMAD4
NCBI Gene ID:	4089
Gene Name:	SMAD family member 4
Gene Aliases:	DPC4, JIP, MADH4, MYHRS
Chromosome Location:	Chr. 18: 51030213 - 51085042 on Build GRCh38
Species:	Homo sapiens

Interrogated Sequence	Translated Protein	Targeted Exon(s)	siRNA Location
NM_001407041.1	NP_001393970.1	9	1165
NM_005359.6	NP_005350.1	9	1580
NR_176264.1	-	10	1642
NM_001407043.1	NP_001393972.1	9	1580

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siRNA Validation Data	
siRNA ID:	s8403
NCBI Gene ID:	4089
Transcript(s):	NM_001407041.1 NM_005359.6 NR_176264.1 NM_001407043.1 NR_176265.1 NM_001407042.1
Validation Results:	83%(+/-1% mRNA knockdown efficiency)(HeLa cell line)
Validation Procedure:	HeLa cells were reverse transfected with 5nM gene-specific <i>Silencer</i> [®] Select siRNA or 5nM <i>Silencer</i> [®] Select Negative Control #2 siRNA. 48 hours post-transfection, gene expression analysis was performed using the TaqMan [®] Gene Expression Cells-to-CT [™] Kit and appropriate TaqMan [®] Gene Expression Assays. 18S rRNA was used as an endogenous control. Validation results are expressed as average percent mRNA remaining, relative to Negative Control siRNA-treated samples, along with the standard deviation of multiple independent experiments.

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☐	Assay ID	Gene	Transcripts	Species	Transcript Type	Product Type	Specifications	Price (JPY)
☐	s8403	SMAD4	6	Human	Coding	Silencer Select	Desalted 5 nmol	定価価格 Each CAT # 4390824 Estimated ship by: 4-5 business days
★ Validated								
View Details Genomic Map Validation Data Recommended Controls Recommended Products								

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Transfection Reagents		
Product Name	Size	Your Price
Lipofectamine™ RNAiMAX Transfection Reagent	1.5 mL	148,400 Each CAT # 13778150
Neon™ Nxt Electroporation System 100-µL Kit with 1-Channel Tubes	96 x 2 reactions	251,000 Each CAT # N10096
Vivofectamine™ VF232 Liver LNP Composition in Ethanol	1 mL	Price: Contact Us Made to order CAT # VF232LVCE
Opti-MEM™ I Reduced Serum Medium	100 mL	3,400 Each CAT # 31985062

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© 2026 Thermo Fisher Scientific Inc. All rights reserved.

All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

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