

INNOVATOR INSIGHT

Platform optimization for efficient AAV purification: insights from a CDMO

Vincent Ravault & Nicolas Laroudie

Over the last decade, the number of clinical trials involving recombinant adeno-associated viral (AAV) vectors has dramatically increased, the diversity of serotypes has expanded, and the demand for larger quantities of highly purified material manufactured to cGMP standards has rocketed. For contract development and manufacturing organizations (CDMOs) like Yposkesi, the manufacturing challenges are centered around flexibility, robustness, and productivity, especially with regards to purification. Universal tools able to address any serotype with minimal process adjustments are critical. In this article, we describe how POROS™ CaptureSelect™ AAVX resin can be used as a pan-affinity tool for the universal capture of AAV vectors, and how Yposkesi optimized the operational parameters to make the resin an efficient, robust, and productive purification platform that fits within the constraints faced by CDMOs.

Cell & Gene Therapy Insights 2022; 8(1), 1–14

DOI: [10.18609/cgti.2022.001](https://doi.org/10.18609/cgti.2022.001)


As a full-service CDMO for innovative gene therapy products, Yposkesi supports customers from early-stage development, including process and analytical development, through to large-scale production and commercial supply of gene therapy products.

Yposkesi produces recombinant adeno-associated virus (rAAV) and recombinant lentivirus

(rLV) vectors using adherent- and suspension-adapted cell platforms. The manufacturing platform at Yposkesi currently includes four independent production suites equipped with 200 L single-use bioreactors, which will evolve to include a 1000 L single-use bioreactor from 2023. Yposkesi is currently building an additional 5,000 m² clinical/commercial

► FIGURE 1

AAV manufacturing process at Yposkesi.

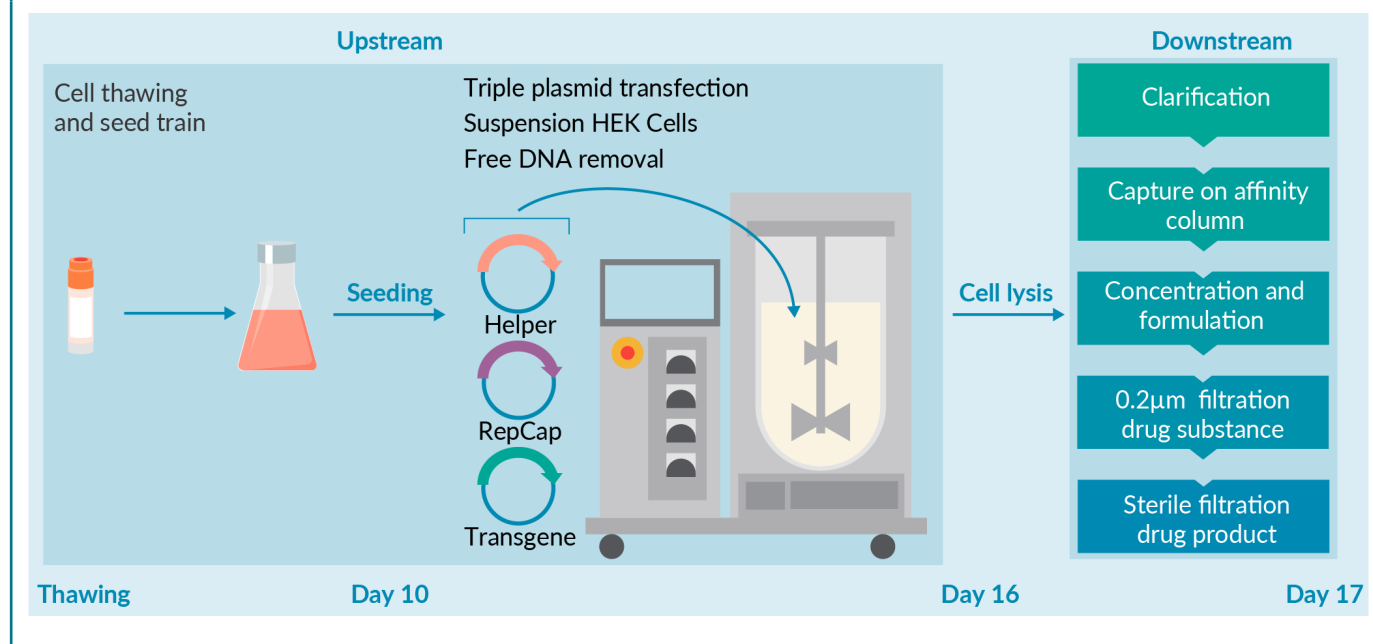

manufacturing plant to support the growing demand for viral vector supply.

This article describes how Yposkesi developed an AAV purification platform for a range of serotypes based on Thermo Fisher Scientific's POROS CaptureSelect AAVX Affinity Resin.

YPOSKESI'S AAV MANUFACTURING PROCESS

The established AAV manufacturing process at Yposkesi is shown in **Figure 1**. AAV vectors are produced by triple plasmid transfection in human embryonic kidney (HEK) cells. The lysate is clarified and then directly loaded onto an affinity column. The eluted vectors are concentrated and formulated, before being sterile filtered. The full process lasts 17 days, from cell thawing to drug product filling.

The current AAV purification process involves the use of different affinity sorbents according to the AAV serotype to be produced. The POROS CaptureSelect AAVX Affinity Resin leans on the use of a ligand derived from a heavy-chain antibody that can bind AAV serotypes 1–9 and synthetic or recombinant AAV vectors, offering a great opportunity to

develop the next AAV purification platform at Yposkesi (**Figure 2**) [1].

EVALUATION OF DYNAMIC BINDING CAPACITY

As a first step to evaluate the AAVX resin as a platform purification solution, the dynamic binding capacity was evaluated using an AAV8 serotype. The binding capacity was assessed using 1 mL-prepacked columns, packed with either POROS CaptureSelect AAV8 or POROS CaptureSelect AAVX. The binding capacity was assessed at 1 and 3 mins residence time on two different feedstocks, each with different initial virus titers.


Clarified supernatant containing AAV8 vectors was directly loaded on the affinity columns until a 10% breakthrough in AAV8 was observed in the flowthrough.

Multiple fractions (column volumes [CV]) were collected at the outlet of the column during the loading phase, and the quantity of capsids was determined by ELISA assay in each collected fraction. The results for the 3 mins residence time are presented in **Figure 3**.

No breakthrough was observed on the AAVX resin at loading volume of up to 1,500

FIGURE 2

Key features of camelid-derived, recombinant expressed ligands used in CaptureSelect™ Affinity Resins.

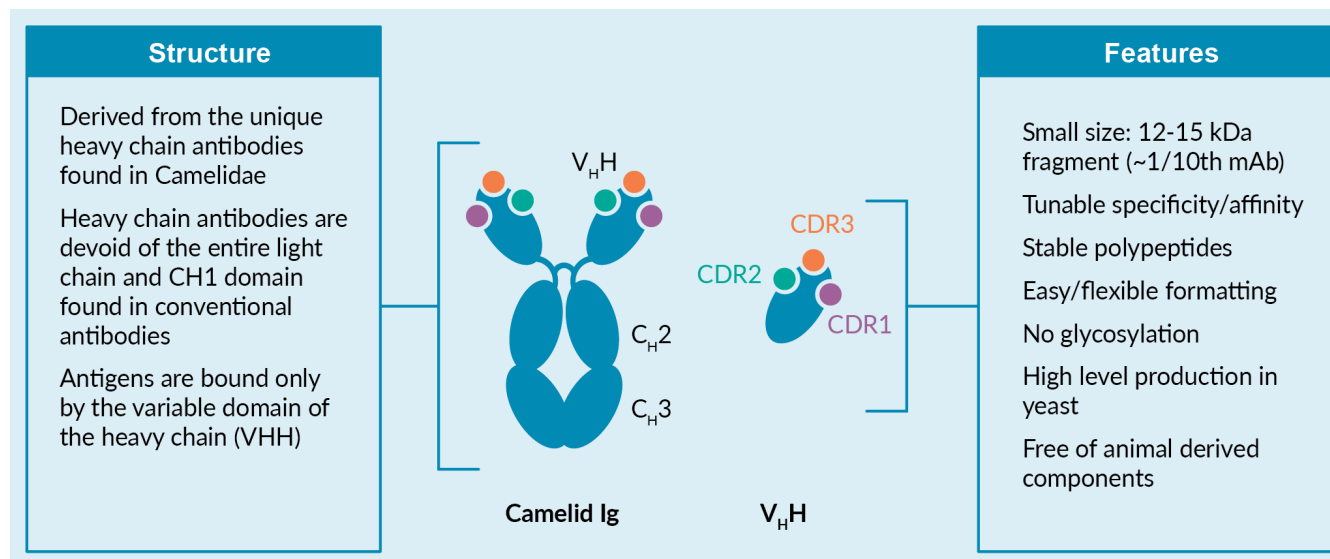


or 2,000 column volumes for the low viral titer and higher titer feedstock, respectively. Both resins showed higher binding capacity when feeds contained a higher vector titer, but overall, the AAVX resin showed a higher

binding capacity for AAV8 than the Poros CaptureSelect AAV8 resin.

Figure 4 shows the binding capacity of AAV8 vectors measured at 1 min residence time on both resins, showing similar binding

FIGURE 3


Binding capacity at 3 mins residence time and 10% breakthrough for POROS CaptureSelect AAV8 and POROS CaptureSelect AAVX at AAV titers of 4.3×10^{10} caps/mL (left) and 4.9×10^{11} caps/mL (right).

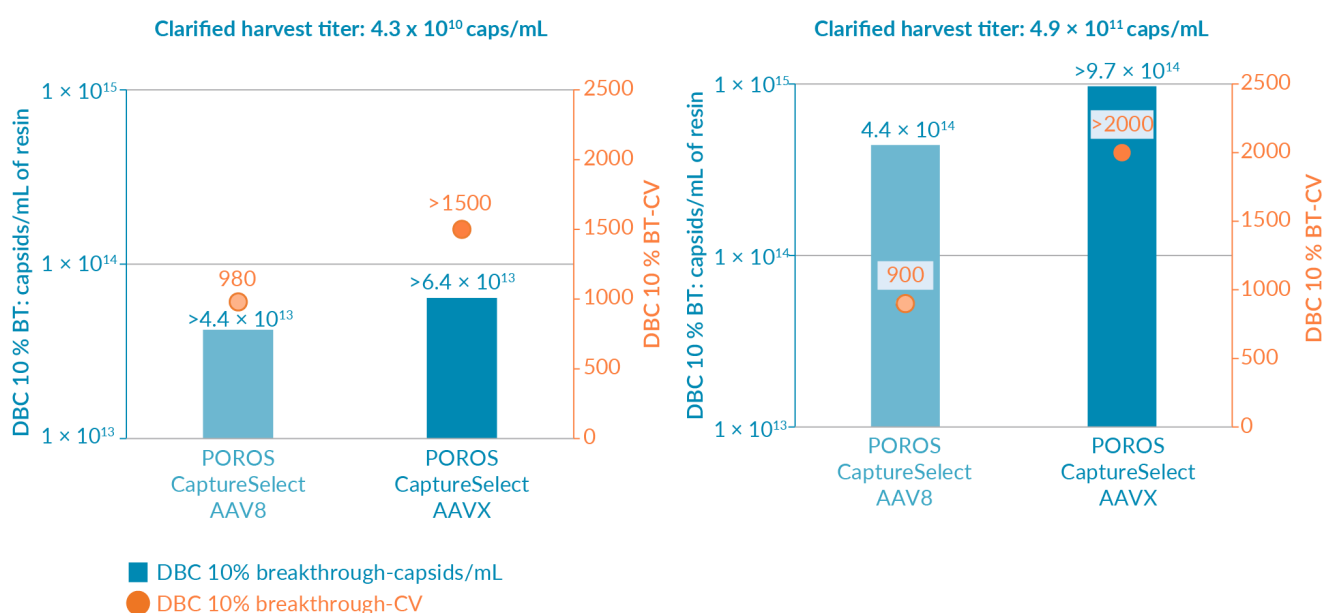

The blue bars represent the binding capacity in terms of capsids per mL of resin at 10% breakthrough. The orange dots represent the column volumes that lead to 10% breakthrough.

► FIGURE 4

Binding capacity at 1 min residence time and 10% breakthrough for POROS CaptureSelect AAV8 and POROS CaptureSelect AAVX at an AAV titer of 5.7×10^{10} caps/mL.

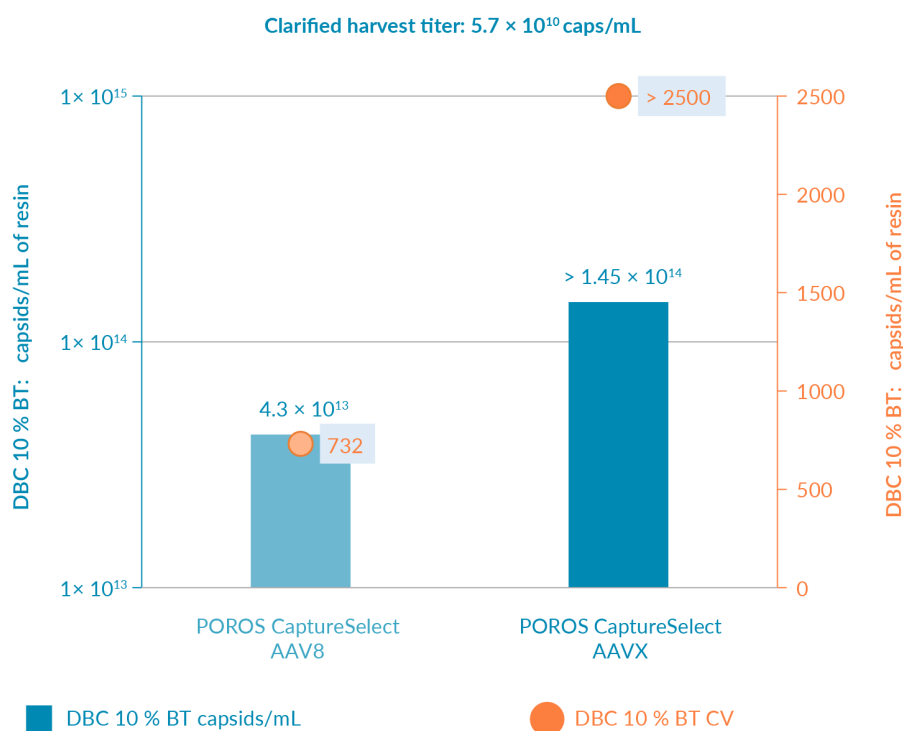

The blue bars represent the binding capacity in terms of capsids per mL of resin at 10% breakthrough. The orange dots represent the column volumes that lead to 10% breakthrough.

capacities compared with the 3 mins residence time. At both residence times, there was no breakthrough on AAVX, with loading volumes up to 2,500 column volumes (CV).

The results from these binding capacity studies led to three main conclusions:

1. The AAVX resin has a better AAV8 binding capacity than the AAV8 resin
2. Binding capacity increases with harvest titer
3. Residence time has no significant effect on the binding capacity

DEFINING OPERATING CONDITIONS FOR PURIFICATION OF AAV8 & AAV2

The operating conditions for the capture of the AAV8 serotype were defined according to

the DBC data obtained previously. Screening of capture conditions was performed on 1 mL-pre-packed columns with AAV8 or AAVX resin (Figure 5).

The material loaded onto the columns was a clarified supernatant containing AAV8 vectors. To align with our AAV manufacturing operating conditions, the maximum loading time selected was 18 hours – (overnight loading). Two residence times were evaluated: 3 mins and 1 min.

The loading volumes selected were 356 CV (for AAV8 and AAVX) with a 3 mins residence time, and 600 CV (AAV8) and 1080 CV (AAVX) with a 1 min residence time. These CVs are all below the resin binding capacities at 10% breakthrough defined earlier (Figures 3 & 4).

After loading and washing, purified product was recovered during the elution step at low pH and was immediately neutralized.

The clarified harvest and eluent were tested for viral genome (VG) titer.

Similar quantities of AAV vectors were loaded on the AAV8 and AAVX resins at 3 mins residence time. As shown in **Figure 6**, the quantity of AAV8 vector recovered after elution and the AAV8 yield was very similar for both resins. The resins showed no significant difference in performance when loading at 3 mins residence time or at 18 hours loading time.

Results at 1 min residence time are shown in **Figure 7**. As a consequence of the different binding capacities at 1 min residence time, the loading times were different for the two sorbents – 10 hours for the AAV8 resin and 18 hours for the AAVX resin. Thus, the total quantity of AAV8 capsid loaded on the resins was around 1.8 times higher for the AAVX resin compared with the AAV8 resin. As a result, the quantity of purified recovered product for AAVX was approximately 1.7 times higher. The step yields for both resins were also very similar and close to 90% which is higher than the yield of around 70% obtained with a residence time of 3 mins.

These results indicate that it will be possible to switch from POROS CaptureSelect AAV8 to AAVX for the purification of AAV8 serotype.

Based on these results with POROS CaptureSelect AAV8, the AAVX resin was also evaluated for the capture of another serotype of AAV: AAV2 (**Figure 8**). The aim was to compare POROS CaptureSelect AAVX with an affinity resin from another supplier, which is currently used at Yposkesi for AAV2 processes.

The screening of the capture conditions was performed on 1 mL prepacked columns. Two residence times were applied for the AAVX resin: 3 mins and 1 min. The residence time applied to the other affinity resin was 8 mins, according to supplier's recommendation.


Three purification conditions were screened for the capture of the AAV2 vector. At 3 mins residence time, the volume of clarified harvest loaded on the column was 455 CV, whereas at 1 min residence time the volume loaded on the column was 1,440 CV. The same starting material was used for all trials.


For the other resin, only 340 CV were loaded since the residence time applied was higher. After column washing, the product was eluted at low pH, and the loading and elution fractions were tested for VG titer.

Using the AAVX resin and decreasing the residence time from 3 min to 1 min resulted in an increase in VG yield from 57% to 89% (**Figure 9**). Using the affinity resin from another supplier with higher residence time

► FIGURE 5

Experimental plan for the definition of operating conditions for purification of AAV8.

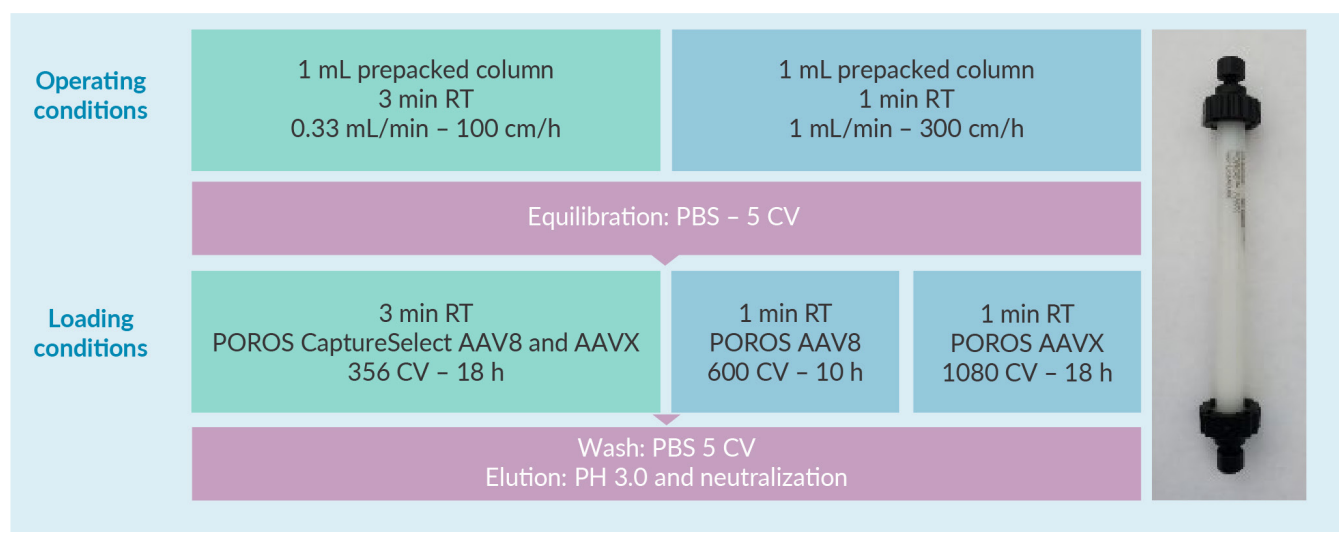

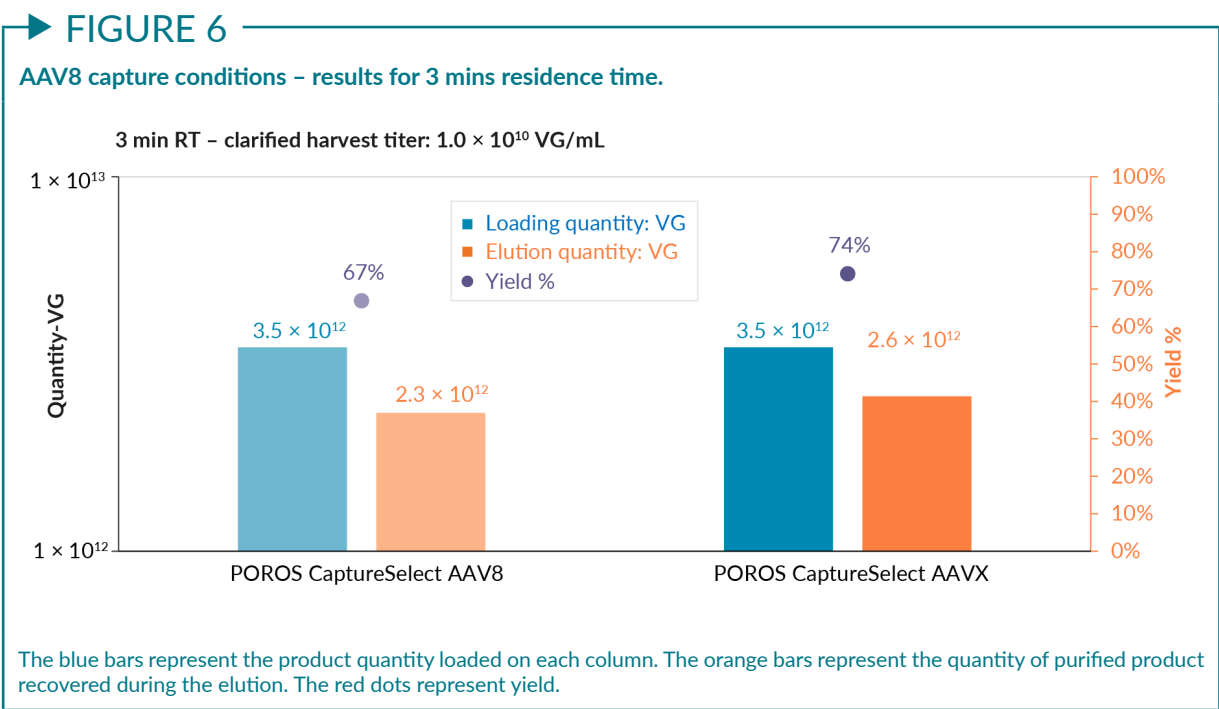


(8 mins, imposed because of the compressibility of the media, and as recommended by the supplier), resulted in a low volume loaded on this column. The AAV2 yield is significantly lower than the yield obtained with AAVX: 48% yield, versus 70–90% yield obtained with AAVX.

This part of the study demonstrated that using a lower residence time results in higher AAV binding capacities for both Thermo Fisher Scientific resins, and that the AAVX


resin shows better results for the capture of AAV8 and AAV2 vectors. The volumes of clarified harvest that can be loaded on AAVX without any AAV breakthrough in the flowthrough are 1080 CV for AAV8, and 1440 CV for AAV2.

The promising results obtained with POROS CaptureSelect AAVX led us to select this resin for the next part of the study and to work with a residence time as close as possible to 1 min.

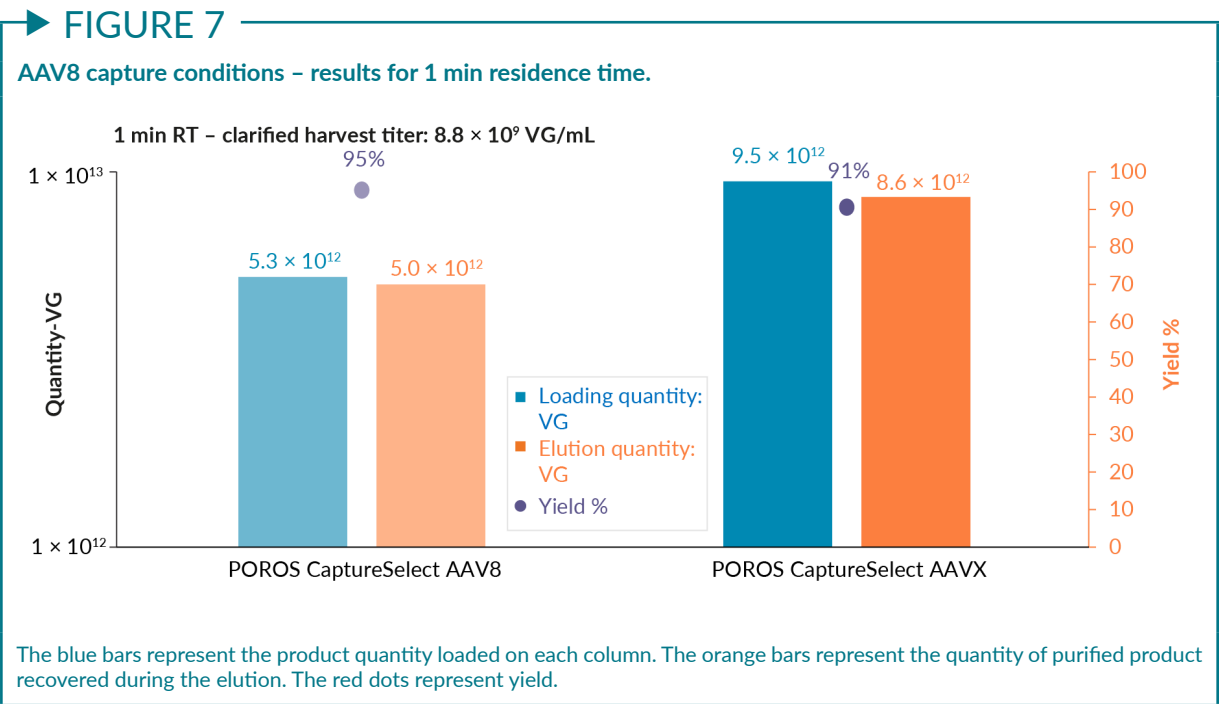

► FIGURE 8

Experimental plan for the definition of operating conditions for purification of AAV2.

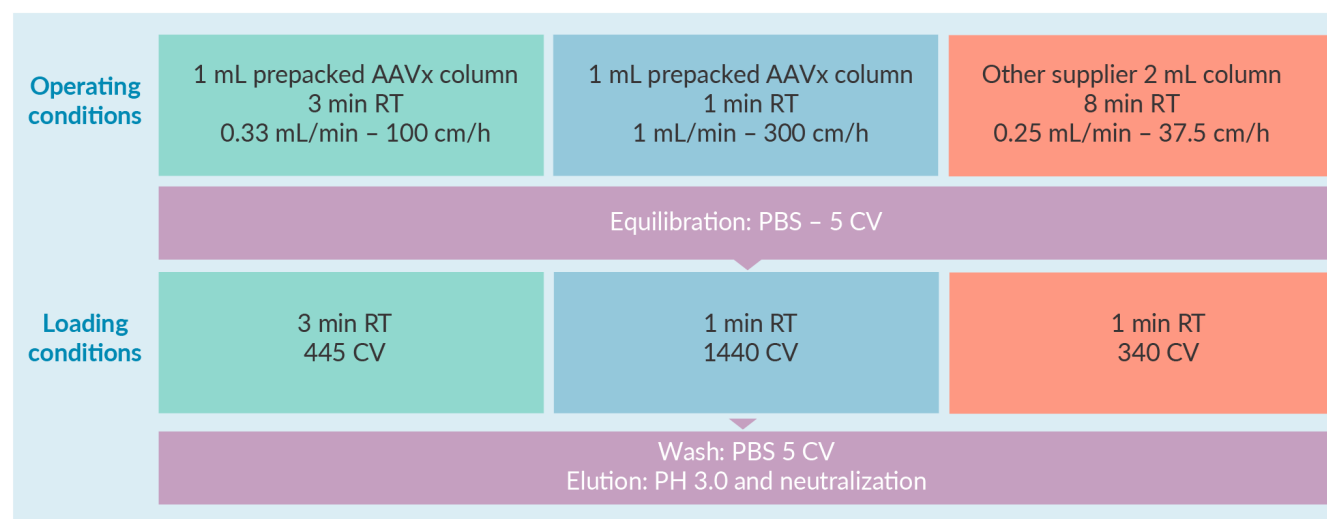


SCALE-UP OF THE CHROMATOGRAPHY STEP

The experimental conditions determined using AAVX for the capture of AAV2 and AAV8 vectors were adapted for the purification of clarified harvest from a 10-liter bioreactor (Figure 10).


The volume of resin necessary to purify a 10 L clarified harvest was calculated by applying the column loading capacity in terms of CV determined previously during the screening for AAV8 and AAV2 processes. This AAVX resin volume was found to be 14.3 mL for AAV8 capture and 11.3 mL for AAV2 purification process.

► FIGURE 9

Definition of the operating conditions for the purification of AAV2 using POROS CaptureSelect AAVX (left) or alternative supplier's affinity resin (right).

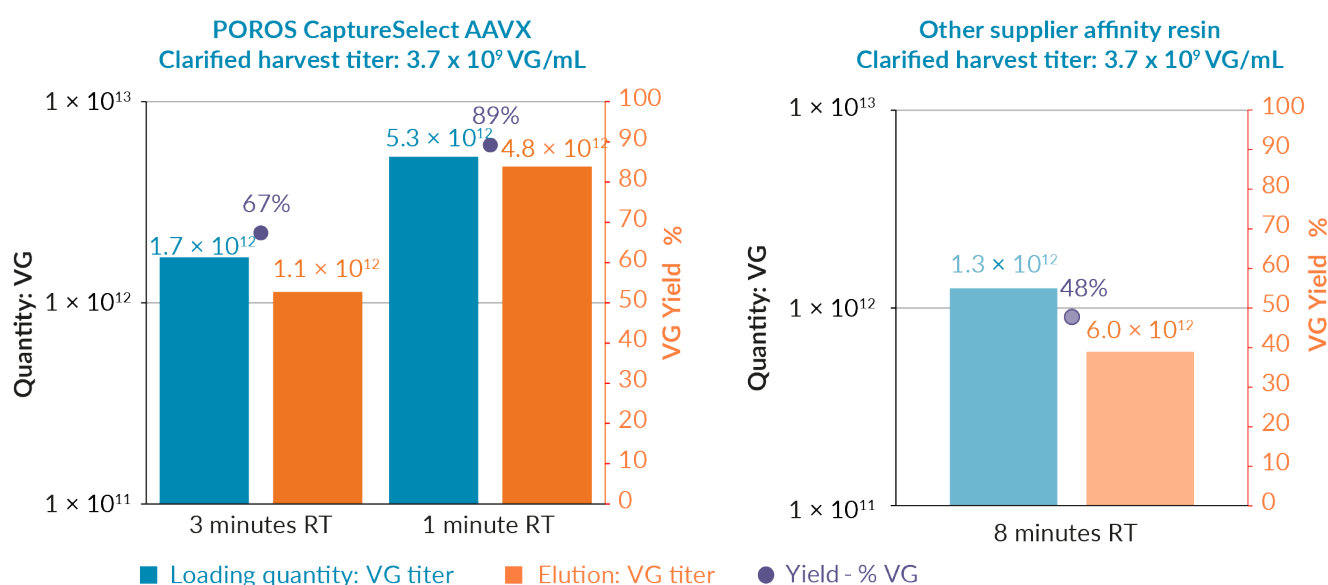

The blue bars represent the product quantity loaded on each column. The orange bars represent the quantity of purified product recovered during the elution. The red dots represent yield.

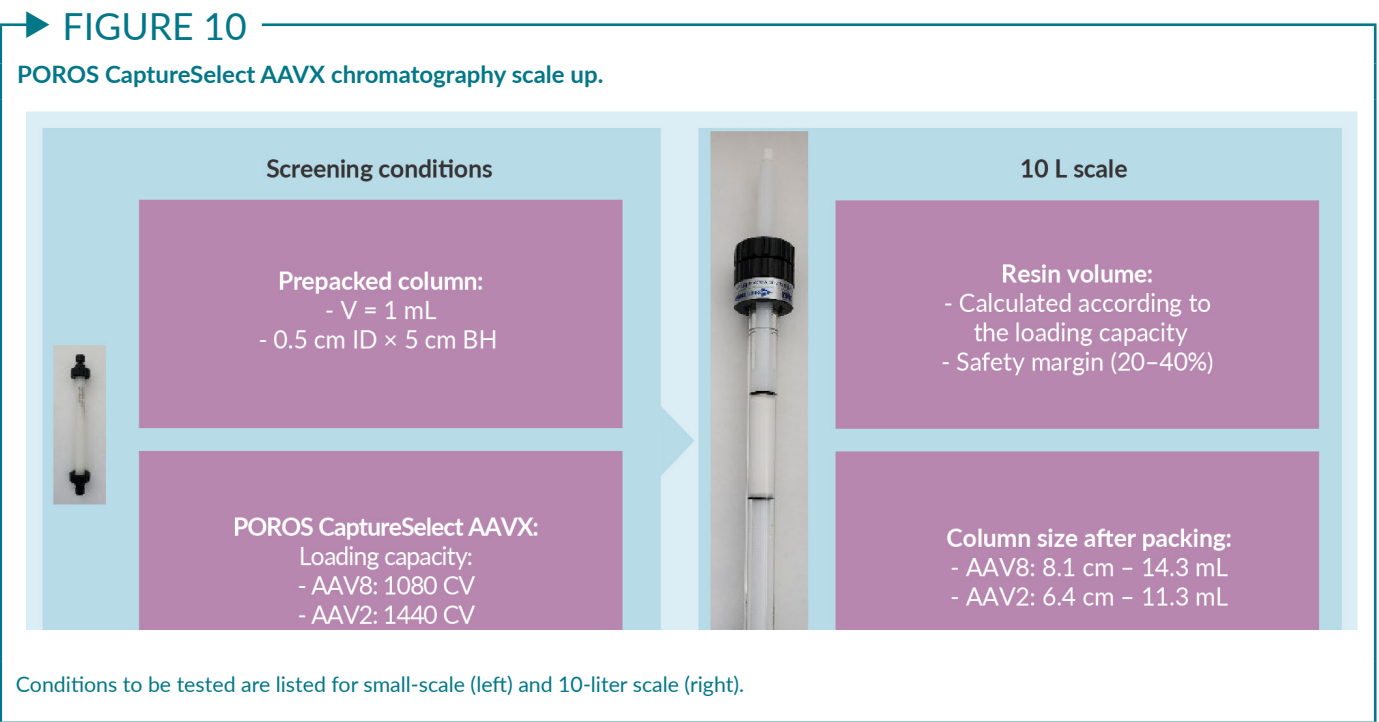

AAVX resin was packed in a 15 mm internal diameter glass column, which allowed for a resin bed height that would be easily transferrable to GMP scale. The column bed height was 8.1 cm for AAV8 purification and 6.4 cm for AAV2 purification.

In order to obtain the starting material for resin evaluation, two 10 L bioreactors were used to produce AAV2 and AAV8 vectors from HEK cells.

After AAV production, cells were lysed, and the lysate was clarified and filtered using a 0.22 µm filter. After lysate filtration, the pool titer was 1.10×10^{11} VG/mL for AAV8 vectors and 3.70×10^9 VG/mL for AAV2 vectors.

The selected operating conditions for the AAVX resin to purify AAV8 and AAV2 from a 10 L clarified harvest are shown in **Figure 11**.


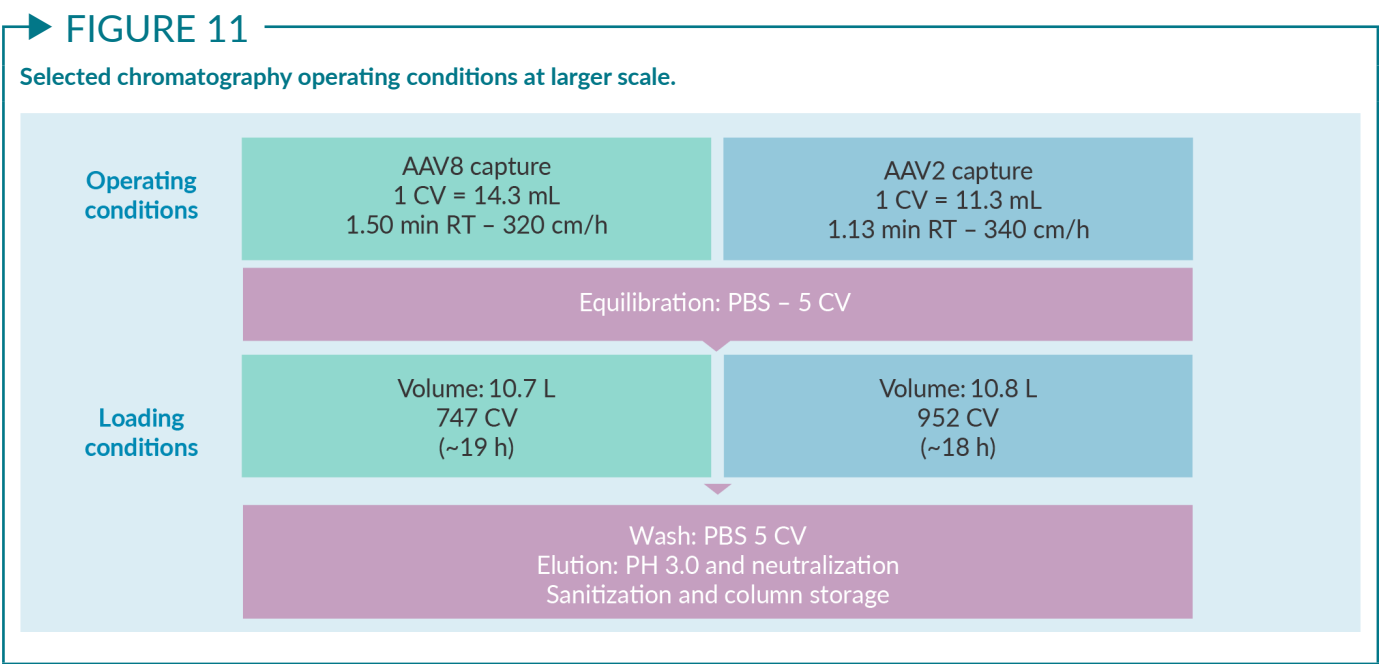


FIGURE 12

AAV capture on POROS CaptureSelect AAVX – yields for AAV8 (left) and AAV2 (right).

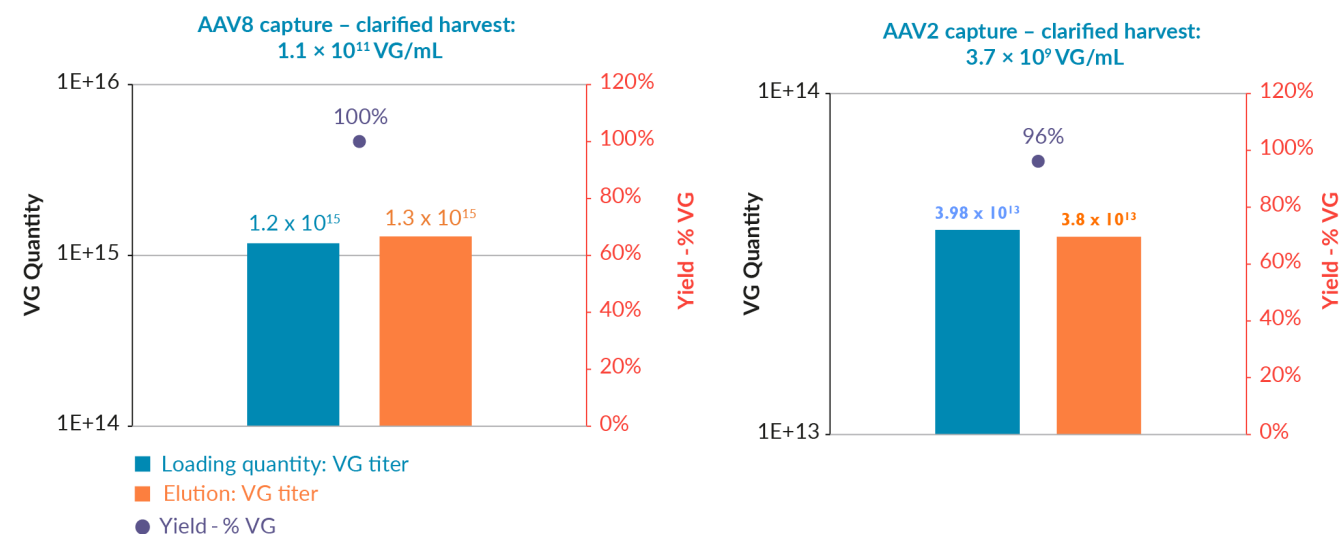

The blue bars represent the product quantity loaded on each column. The orange bars represent the quantity of purified product recovered during the elution. The red dots represent yield.

The residence time for both AAV processes was close to 1 min. 747 CV of clarified harvest were loaded on to the AAVX resin for AAV8, and 952 CV for AAV2, while the loading times were in the same range. The purified products were recovered during elution at low pH and then neutralized. VG titers, total protein content, and residual DNA levels were assayed in the clarified harvests (starting materials) and in the elution fractions.

The pressure was monitored at the inlet of the column during the loading step for the AAV8 and AAV2 capture process. The pressure slightly increased during the loading stage but stayed within an acceptable range. The pressure was around 1.5 bars at the end of the loading step, which helps to provide good conditions for a transfer to GMP scale.

Even though VG titers in the starting material were very different for the AAV2 and AAV8 serotypes, the final yields of the capture step are close to 100% for both serotypes and there was good scalability from lab scale development to the 10 L scale (Figure 12).

Additional experiments revealed that the purity of AAV vectors captured with AAVX

resins appears to be very high. There was an impurity reduction of over 99% in the purified product after capture on AAVX for each serotype. This clearance rate could be even further optimized by adding an intermediate washing step or implementing a polishing column after the AAV capture step.

CONCLUSIONS

This long-term study with POROS CaptureSelect AAVX resin has highlighted several advantages of AAV capture using this resin compared to other affinity resins commercially available:

- ▶ Flexibility in terms of serotypes: capture of AAV1 to AAV9 serotypes and synthetic and recombinant serotypes
- ▶ Possibility to standardize a purification platform for several AAV serotypes with only a few adjustments
- ▶ Cost reduction due to shorter residence times and very high loading volumes.

- ▶ Low level of impurities captured on the resin. This could be further optimized for each serotype if needed (wash conditions screening or addition of a polishing step)
- ▶ Good scalability of the downstream platform. It is compliant for a large-scale GMP AAV manufacturing process

Overall, Yposkesi concluded that the POROS CaptureSelect AAVX resin appears to be a great tool to improve purification processes in terms of quality, cost, and standardization. Yposkesi plan to implement this resin for the purification of other AAV serotypes.

ASK THE EXPERTS



Nicolas Laroudie (Thermo Fisher Scientific) joins **Vincent Ravault** (Yposkesi) to answer readers' questions on implementing POROS CaptureSelect technology into viral vector production.

Q Can the POROS CaptureSelect resin be cleaned and re-used?

NL: Yes, the resin can be cleaned and reused. Many customers use the resin once, particularly CDMOs that deal with multiple serotypes and multiple transgenes and want to avoid spending a lot of time validating cleaning. But the resin can absolutely be cleaned and reused, and many customers are doing that.

Notably, the resin is not very caustic stable, and so for cleaning, we do recommend using acidic solutions such as phosphoric or citric acid. In case of very dirty resin, we advise additional cleaning with chaotropic agents, such as guanidine hydrochloride or urea.

I would encourage people who want to clean and re-use the resin to reach out to their local application specialist, who can help them develop a process for this.

Q Apart from AAV2 and AAV8, do you have experience with other AAV serotypes and POROS CaptureSelect AAVX?

VR: The goal for Yposkesi now is to expand this platform to a broad range of AAV serotypes. With our experience of AAV2 and AAV8 serotypes, we know that the AAVX resin is a good solution to use as a purification platform.

Currently, we are working with AAV5, 6, and 9, and the results so far are promising. We also know that we can work with modified capsids.

Q Is the resin available in a pre-packed format?

NL: Yes, we do have pre-packed formats. We have 1 and 5 mL pre-packed formats available that are compliant with standard benchtop chromatographic systems. We also have robocolumns available, at 200 µL and 600 µL, for high-throughput screening.

Of course, the resin can be purchased as bulk material and our local Field Application Specialists are happy to support customers in packing the resins in their own columns, whatever the scale.

Q Which additional washing conditions would be suitable for host cell protein and host cell DNA reduction?

VR: Several washing conditions are interesting to assess. For example, you can add an extra washing step using high salt concentration. You can also wash your column with a low pH buffer in order to remove impurities from the column before recovery of AAV in the eluate.

If you decide to implement the second washing step you have to be careful that your washing condition won't affect the integrity of your capsid. Moreover, if the washing step pH is too close to the pH of the elution buffer, a significant quantity of capsids could be eluted during your washing step, and as a result, lower your AAV yield during the elution.

Q Is the resin GMP compliant?

NL: The resin is used in GMP manufacturing by many of our customers. While not manufactured in a cGMP process, the resin is produced under an ISO 13485 environment.

When you purchase the resin, you can request the regulatory support package, including documents regarding quality, stability, production, control method, and so on. Those documents are useful when you make a product and submit a dossier to a regulatory agency.

For each of our commercially available CaptureSelect resins, we developed an ELISA assay to monitor the level of ligand leakage over the purification process.

Q Vincent, why did you use two different analytical methods during this study – ELISA for dynamic binding capacity and viral genome titers at termination during your screening?

VR: During our DBC study, a lot of fractions were collected in the flowthrough at the outlet of the column in order to calculate 10% breakthrough for AAV vectors, so we needed to use a high throughput assay for the analysis of the first full fractions. The ELISA assay allowed us to test several samples in parallel and to get the results quickly, in around half a day.

In the screening study, the number of samples was much lower – only two samples for each set of conditions screened were produced – so here we used an internal assay for the quantification of the viral genome titer in the product. The viral genome titer was determined by qPCR for each serotype.

Q Which resin can be implemented for a polishing step?

VR: Several different resins can be implemented for this step. Commonly, an anion exchanger is implemented in order to reduce host cell protein and host cell DNA. Anion exchange also has the capability to separate empty and full AAV capsids, and some suppliers have developed resins specifically for the polishing step. For more information, you can contact chromatography resin suppliers.

REFERENCE

1. Sierkstra L. Overcoming purification challenges in antibody therapeutics manufacturing with subdomain-specific affinity solutions. *Immuno-Oncology Insights* 2021; 2: 229–39.

BIOGRAPHIES

Vincent Ravault

DSP Expert, Process Development and Industrialization Department, Yposkesi

Vincent Ravault is a technician within the Process Development and Industrialization department at YposKesi since 2018. He focuses on upstream and downstream process development and optimization for AAVs and Lentiviruses vectors to support and advance their large-scale production for gene therapies. He is also involved in the technical transfer of processes for viral vectors from pilot scale to manufacturing scale. Graduated of a Biochemistry degree, Vincent Ravault has over 15 years of experience at Pall Life Sciences where he was in charge of the technical support for chromatography resins. His main role at Pall was to provide purification solutions and strategies to customers. He was also involved in the promotion and evaluation of new products for downstream.

Nicolas Laroudie

Staff Scientist, Field Applications, Thermo Fisher Scientific

Biochemist by education, Nicolas Laroudie used to work for Généthron, France, between 2001 and 2011 as Head of Downstream Development. He was leading a team in charge of developing and scaling-up purification processes for AAV, retroviral and lentiviral vectors

used in gene therapy treatments. He then joined Merck Millipore as a BioManufacturing Engineer where he used to technically support European customers for all DSP technologies - from clarification to sterile filtration, including TFF and systems - with a strong focus on chromatography. In particular, he took an active role in the establishment of a fully continuous, large-scale disposable DSP process for the purification of a monoclonal antibody, within the framework of a large multi-company European consortium. He eventually joined ThermoFisher Scientific in 2019 as Field Application Specialist for purification, technically supporting the implementation of POROS and CaptureSelect chromatography products for south-western European customers.

AUTHORSHIP & CONFLICT OF INTEREST

Contributions: All named authors take responsibility for the integrity of the work as a whole, and have given their approval for this version to be published.

Acknowledgements: None.

Disclosure and potential conflicts of interest: The authors declare that they have no conflicts of interest.

Funding declaration: The authors received no financial support for the research, authorship and/or publication of this article.

ARTICLE & COPYRIGHT INFORMATION

Copyright: Published by Cell and Gene Therapy Insights under Creative Commons License Deed CC BY NC ND 4.0 which allows anyone to copy, distribute, and transmit the article provided it is properly attributed in the manner specified below. No commercial use without permission.

Attribution: Copyright © 2022 Thermo Fisher Published by Cell and Gene Therapy Insights under Creative Commons License Deed CC BY NC ND 4.0.

Article source: This article is a transcript of a webinar, which can be found [here](#).

Webinar recorded: Oct 26 2021; **Revised manuscript received:** Jan 4 2021; **Publication date:** Jan 24 2022.


ThermoFisher
S C I E N T I F I C



This is a transcript of a webinar.
You can also watch the recorded webinar:

WATCH NOW



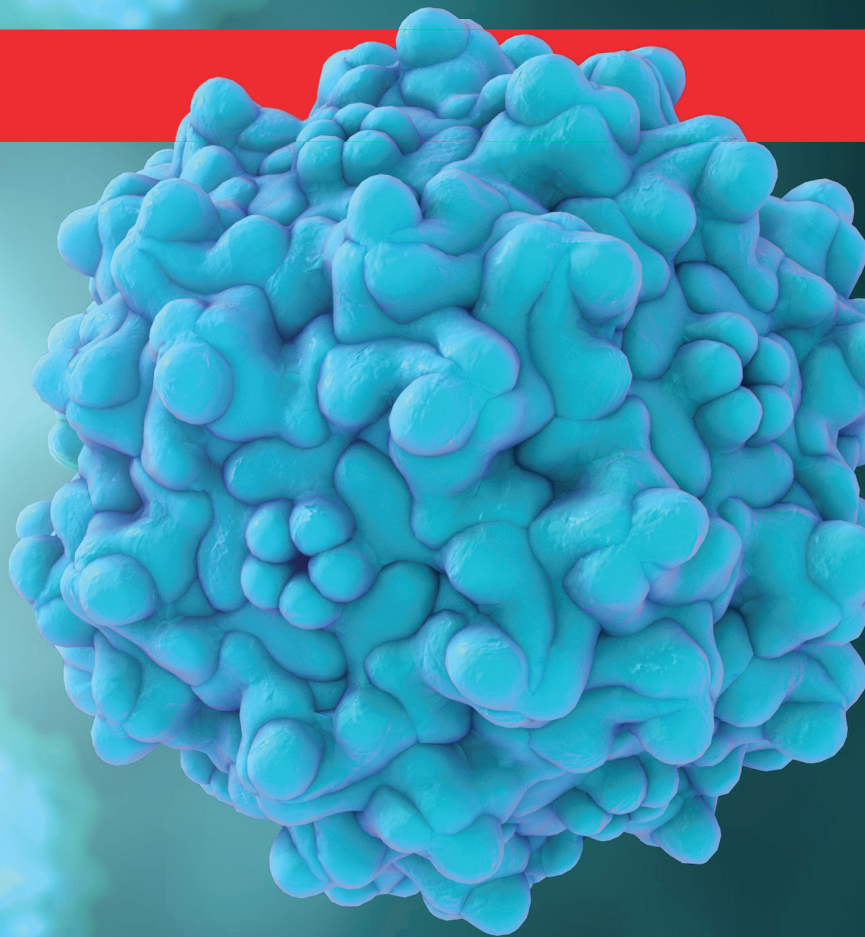

Accelerate advancement in gene therapy

Improve productivity of your AAV downstream process

Reduce purification steps and increase throughput in your viral vector purification workflow. Thermo Scientific™ POROS™ CaptureSelect™ AAV affinity resins allow you to:

- Maximize downstream process efficiency
- Achieve high purity and yield in a single purification step
- Scale up to clinical and commercial gene therapy manufacturing



To learn more, visit thermofisher.com/aav-purification

ThermoFisher
SCIENTIFIC

For Research Use or Further Manufacturing. Not for use in diagnostic procedures. © 2020 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. COL23924 0220