


WildFire: A Simple Monoclonal Colony Generation Technology without Emulsion PCR

Zhaochun Ma, Raymond Lee, Jonathan Erikson, Swati Goyal and Kai Lao,
Life Technologies, 850 Lincoln Centre Drive, Foster City, CA 94404, USA

ABSTRACT / INTRODUCTION

We have developed a new massive parallel monoclonal template generation technology, called "WildFire", where sequencing libraries are in-situ isothermally amplified directly on the surface of a 5500 flowchip. Sequencing libraries are added directly to the 5500-series Genetic Analyzer flowchip, whose surfaces have been coated with a special library-adaptor capture oligonucleotide. A DNA polymerase reaction mix is added, and in a single isothermal step for 30 minutes, single templates are amplified in-situ on the flowchip. The net density of sequencing-colonies created in this manner far exceeds anything currently utilized in next-generation sequencing, reaching 1.1 million colonies per mm² per flowchip surface. During in-situ amplification, the capture oligonucleotide is "consumed", and each individual nucleic-acid fragment "spreads" (like a WildFire) inside the flowchip until reaching an adjacent library fragment "meet", the amplification step terminates, because all of the surface-bound primer was consumed. These "self-assembled", spatially resolved, monoclonal colonies, are then sequenced by SOLiD™ chemistry. The resulting colony-sequencing reads maintain the same high accuracy as our bead-based method. Full genomes (from bacterial to human), exomes (human), and transcriptomes (human) have now been sequenced using WildFire technology. Here we also demonstrate a novel reverse sequencing method with only SOLiD™ forward probes, called WildFire Paired-End sequencing (wPE) using a template switch strategy.

Radically Simplified 5500 Sequencing Workflow

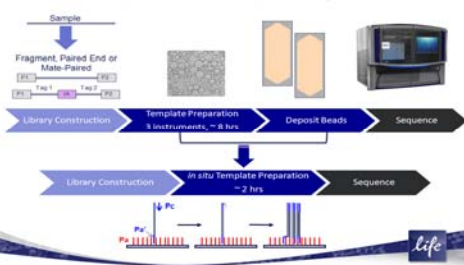

MATERIALS AND METHODS

All targeted capture enrichments (singleplex and multiplex) were done with Ion TargetSeq™ Exome Enrichment Kit probe-sets according to the manufacturer's protocols. DNA was obtained from the Coriell Institute and 5500xl W, 5500xl Genetic Analyzer libraries were constructed using kits from Life Technologies using 5500 SOLiD® Fragment Library Barcode Adaptors 1-16 Kit with and without the 5500xl W conversion module respectively. Two types of library samples including NA12878 and HuRef were used in the preparation of 4 barcoded libraries using an AB Library Builder™ instrument, followed by pippin prep for size selection. This was followed by the Ion TargetSeq™ (v2) exome enrichment. An aliquot of the library pool was used for template preparation on FlowChip using the Wildfire "template walking" protocol. Sequencing on 5500xl W Genetic analyzer was done in single-lane runs for 50x50 paired-end tag. Data analysis, including mapping and targeted resequencing analysis and variant calling, was done with LifeScope™ software.

RESULTS

Figure 1. 5500xl W in-situ Template Preparation Is a Fast Isothermal Amplification Generating Ultra High Density Colonies

Isothermal Template Walking (TW)



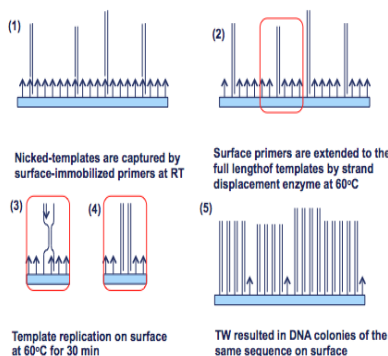


Figure 2. 5500xl W Colonies "Self-Assemble" on Both Sides of Flow-Chip, Further Doubling Output and Reducing the Reagent Cost

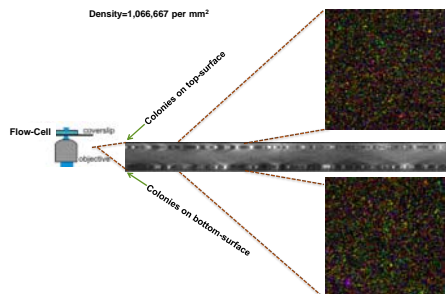

Note: dual surface format available outside US only

Figure 3. WildFire Pair-end Reads(wPE) by template switch

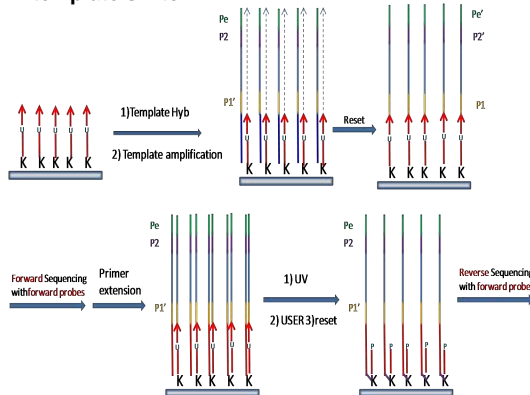

K: photo cross-linking reagent, U: deoxyuridine

Table 1. A Wildfire 50x50 Paired End (wPE) Run Generated 21.3 Gb Mapped /Lane (570 top only layer panels)

Filtered Throughput	38 GB
Mapped Throughput	21.34 GB
Total F3 Tag Mapped	276M
Total F5 Tag Mapped	199M
Throughput (Gb/day)	9.5
Both F3 and F5 Tag Mapped %	55%
F3 Tag Mapped only	73%
F5 Tag Mapped only	53%
Panels scanned	570
Reads in Target	249M (64.92%)
Reads off Target	134M (35.07%)
Percent of target bp not covered	4.3%
Average Depth of Coverage within targets	216X
Percent of target bp covered at >=20X	87.81%
dbSNP heterozygous concordance	96.18%
dbSNP homozygous concordance	99.40%

CONCLUSIONS

WildFire technology greatly improves NGS workflow, increases throughput, and significantly decreases net cost-per-genome. A novel pair-end sequencing method was also developed by using a template switch strategy. With top layer only, a single lane of 50x50bp PE run, we generated over 21Gb mapped data with an average depth of 216X, >95% of target bases covered and >87% covered at 20X depth with a db heterozygous SNP concordance of 96%.

Acknowledgement

We thank Devin Do and Kamini Varma for providing TargetSeq™ exome library.

For Research Use Only. Not intended for any animal or human therapeutic or diagnostic use.

© 2012 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation and/or its affiliate(s) or their respective owners.